Instructions for use

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To Open BioBall® Aseptically remove stopper.

Agar Plate

1. Tip BioBall into the centre of the plate.
2. Rehydrate by pipetting 100 µL of sterile water or 0.9% saline solution directly onto BioBall.
3. Wait 30 seconds for BioBall to dissolve.
4. Use a sterile plastic spreader to evenly spread the dissolved BioBall over the plate surface.
5. Ensure plate is dry before inverting and incubating.

Membrane Filtration

1. Tip BioBall into liquid sample.
2. Mix by repeatedly inverting sample
3. Filter the sample using standard test method.
4. Follow standard test method for incubation

Pour Plate

1. Pipette 1ml of sterile water or 0.9% saline solution into pour plate.
2. Tip BioBall into the sterile water/saline solution.
3. Wait 30 seconds for BioBall to dissolve.
4. Add molten agar and mix according to standard test method.
5. Incubate as required.

Petrifilm®

1. Tip BioBall onto the centre of Petrifilm®.
2. Rehydrate by pipetting 100 µL of sterile water or 0.9% saline solution directly onto BioBall.
3. Follow standard test method.
4. Incubate as required.
5. Petrifilm® is a registered trademark belonging to 3M™ Corporation.

Adding to Samples/Broths

1. Tip BioBall® into sample/broth
2. Mix
3. Incubate as required

Intended Use

BioBall SingleShot is a microbiological reference material containing a precise number of viable bacterial cells. It is designed for use as a quantitative quality control sample.

Description

BioBall SingleShot is a freeze dried water soluble ball containing a precise number of micro-organisms. The ball is a white sphere approximately 3mm in diameter.

Quality Control

BioBall SingleShot delivers a level of precision previously unobtainable in microbiology. Each batch of BioBall SingleShot is quality controlled by testing a minimum of 10% of the batch up to a maximum of 50. Vials are selected at random from throughout the batch and plated onto non-selective agar plates and DRBC for

Aspergillus species. The plates are incubated then counted. Each standard batch of BioBall meets the following criteria (unless otherwise stated):

Mean	28-33 cfu
Standard Deviation	3 cfu or less

Certificate of Analysis

Certificates of Analysis can be downloaded from the BTF website: www.bioball.com. The certificate states the mean cfu and standard deviation of the batch. It also states the 95% Prediction Interval. The standard deviation is a measure of variability within the batch, the mean is the average cfu count tested on nonselective agar and DRBC for *Aspergillus* species. The Prediction Interval is a statistical analysis of the quality data stating that 95% of your expected results will fall within this range. Our team of technical experts can be contacted regarding the interpretation of results. The micro-organisms used in BioBall® SingleShot are derived from known reference strains. Their identity is confirmed by genetic typing. ATCC is a trademark of the American Type Culture Collection.

Limitations

CFU count (mean, CV and standard deviation) stated in this document and BioBall promotional literature, and batch-specific Quantification Data reported on BioBall Certificates of Analysis, are determined on non-selective culture media (unless otherwise stated). Selective media vary greatly in their inhibitory properties amongst different formulations and amongst brands with similar formulations. Recovery rates from BioBall may be reduced relative to the selectivity of the media. Users of BioBall with selective media are advised to establish their own expectations for recovery rate for each BioBall strain on selective media and to base their performance expectations on data derived from the specific formulations and brands of selective media used in their laboratories. Where recovery rates are reduced on very selective media, users should consider using a greater inoculum to meaningfully assess the limits of detection of their media.

Accreditation

BTF as a Reference Materials Producer is the first company worldwide accredited for quantitative microbiological reference standards, including BioBall.

Technical Assistance

BioBall SingleShot can be used for other applications. For technical assistance please contact your bioMerieux or BTF customer support team.

Safety Precautions

BioBall SingleShot contains viable and potentially pathogenic bacteria and should only be handled by experienced laboratory personnel trained in the safe handling of these micro-organisms. The number of micro-organisms contained within BioBall SingleShot is low. However, BioBall SingleShot should be handled as if potentially infectious. Refer to your national safety guidelines. After use, dispose of packaging in accordance with appropriate biohazard disposal practices. Do not use the product if the packaging is damaged. Do not use the product after the expiry date. The product should be used according to the procedure described in this instruction for use document. Any modifications may affect the results.

Warranty

BTF warrants to the original purchaser only that the products will conform to the quantity and contents stated on the product labels and quality certificate for the duration of the stated shelf life.

Limits of Liability

BTF's sole obligation and the purchaser's exclusive remedy under the above warranty is limited either to replacement, at BTF's expense, any products which are defective in manufacture (provided that the purchaser must return the products, transportation prepaid, to BTF) or, at BTF's option, refund the purchase price. All claims must be notified to BTF within 14 days of using the product. The above warranty does not apply to any products which have been altered outside BTF, nor to any products which have been subjected to misuse or mishandling, or use, storage or handling other than in accordance with instructions supplied by BTF. ALL OTHER WARRANTIES, EXPRESS, IMPLIED OR STATUTORY, ARE HEREBY SPECIFICALLY EXCLUDED, INCLUDING BUT NOT LIMITED TO WARRANTIES OR MERCHANTABILITY OR FITNESS FOR A PARTICULAR PURPOSE AND NONINFRINGEMENT. BTF's maximum liability is limited in all events to the price of the products sold by BTF. IN NO EVENT SHALL BTF BE LIABLE FOR ANY SPECIAL, INCIDENTAL OR CONSEQUENTIAL DAMAGES OR LOSS OF PROFITS. The limitations set out above apply to the fullest extent permitted by law. If any provision of these terms and conditions is held to be invalid, illegal or unenforceable, the enforceability of the remaining provisions will be unaffected.

Use and Indemnity

The purchaser assumes all risk and responsibility in connection with the use, storage, handling and disposal of the products. The purchaser must ensure that all use, storage, handling and disposal of the products is in compliance with all applicable laws, regulations and guidelines and in accordance with any instructions supplied by BTF, and must take all appropriate safety precautions. The purchaser indemnifies BTF and its directors, officers and employees and agents from and against all claims, actions, liabilities, costs and expenses arising directly or indirectly from the use, storage, handling and disposal of the products, except to the extent arising directly from the negligence of BTF.

© Copyright 2009 BTF Pty Ltd Sydney Australia exercising jurisdiction there. BioBall[®] does not contain organisms sourced from The American Type Culture Collection[®]. BioMérieux, the blue logo, BTF and BioBall are used, pending and/or registered trademarks belonging to bioMérieux S. A., or one of its subsidiaries. Any other name or trademark is the property of its respective owner.

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© Copyright 2009 BTF Pty Ltd Sydney Australia exercising jurisdiction there. BioBall® does not contain organisms sourced from The American Type Culture Collection®. bioMérieux, the blue logo, BTF and BioBall are used, pending and/or registered trademarks belonging to bioMérieux S. A., or one of its subsidiaries. Any other name or trademark is the property of its respective owner.

Manufactured by:

BTF Pty Ltd.

Sydney Australia

Tel. 61 (0) 2 8877 9150

Fax. 61 (0) 2 8221 9589 www.bioball.com

Sales & General Enquiries:

bioMérieux S.A

69280 Marcy l'Etoile - France

Tel. 33 (0) 4 78 87 20 00

Fax. 33 (0) 4 78 87 20 90

www.biomerieux.com www.biomerieux-industry.com

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© Copyright 2009 BTF Pty Ltd Sydney, Australie, ayant compétence dans ce pays. BioBall® ne contient pas d'organismes issus de l'American Type Culture Collection®. BioMérieux, le logo bleu, BTF et BioBall sont des marques utilisées, déposées et/ou enregistrées appartenant à bioMérieux S.A. ou une de ses filiales. Toute autre appellation ou marque déposée appartient à son propriétaire respectif.
© Copyright 2009 **BTF Pty Ltd** オーストラリア、シドニーが管轄権を有するものです。
© Copyright 2009 BTF Pty Ltd Sydney Australia sob esta jurisdição. A BioBall® não contém microrganismos provenientes da American Type Culture Collection®. A bioMérieux, o logotipo azul, BTF e BioBall são utilizados, pendentes e/ou registados, propriedade exclusiva da bioMérieux S. A., ou de uma das suas filiais. Qualquer outro nome ou marca registada é da propriedade do respectivo proprietário.
© Copyright 2009 BTF Pty Ltd Sydney Australia donde ejerce la jurisdicción. BioBall® no contiene organismos provenientes de The American Type Culture Collection®. bioMérieux, el logo azul y BioBall son marcas comerciales utilizadas, depositadas y/o registradas pertenecientes a bioMérieux S.A. o a alguna de sus filiales. Cualquier otro nombre o marca comercial es propiedad de su dueño respectivo.
© Copyright 2009 BTF Pty Ltd Sydney Australia. BioBall® nie zawiera organizmów pochodzących z American Type Culture Collection®. bioMérieux, jego niebieski znak, BTF i BioBall są używanymi, i/lub podczas rejestracji znakami towarowymi należącymi do bioMérieux S. A. lub jednego z jej przedstawicieli. Jakiegokolwiek inne nazwy i znaki towarowe należą do odpowiednich właścicieli.
© Copyright 2009 BTF Pty Ltd Sydney Australia, joghatóság gyakorlása ott. A BioBall® nem tartalmaz American Type Culture Collection®-ből származó mikroorganizmusokat. A bioMérieux, a kék logó, a BTF és a BioBall a bioMérieux S. A. vagy valamely leányvállalata által használt, bejegyzésre váró és/vagy bejegyzett védjegy. Minden más név vagy védjegy saját tulajdonosának képezi tulajdonát.